

Obesity Hormone/Adipokine ELISA Combo Kit

Catalog # EA-7084

(For Research Use Only)

Introduction

The Signosis Obesity Hormone/Adipokine ELISA Combo Kit offers multiplexed quantification of six major obesity-related hormones and two cytokines—insulin, amylin, glucagon, GLP-1, leptin, adiponectin, IL-6, and TNF α —in a single 96-well plate (12 wells per target). Compatible with serum, cell, tissue lysates, and various biological fluids from human, mouse, and rat samples, this assay enables sensitive, specific, and high-throughput measurement for studies in endocrinology, metabolic disease, and obesity research. The streamlined protocol is ideal for investigating metabolic pathways in diabetes, obesity, and related disorders, providing a comprehensive solution for simultaneous analysis of key metabolic biomarkers.

Principle

Competitive ELISA

The hormones insulin, amylin, glucagon, and GLP-1, are analyzed using a competitive ELISA assay. In this assay, the plate is coated with the target hormone conjugated to a protein, which competes with the free hormones in the samples that are added to the plate. After the samples containing free hormone are added to the plate, an antihormone antibody labeled with biotin is added to the samples and binds to both the free hormones and conjugated hormones immobilized on the plate. The plate is then washed to remove any sample and antibodies not binding to the hormones bound on the plate. The antibodies that are bound to the immobilized hormones on the plate are measured by adding Streptavidin-HRP, which binds to the biotin-labeled antibodies. To detect the HRP, TMB substrate is added, which produces a blue color when it interacts with HRP. After the blue color fully develops, the reaction is terminated with an acidic Stop solution, which changes the blue color to yellow. The intensity of the yellow color is measured at an absorbance of 450 nm and is used to determine the hormone level in the sample. Because this is a competitive ELISA assay, the signal is inversely proportional to the hormone concentration, since higher free hormone concentrations will prevent the labeled antibody from binding to the plate.

Sandwich ELISA

The peptides leptin, adiponectin, IL-6, and $TNF\alpha$, are analyzed using a sandwich ELISA assay. In this assay, peptides present in the sample bind to specific capture antibodies pre-coated on a plate. After unbound components are washed away, biotin-labeled detection antibodies are added to bind the immobilized hormones. Streptavidin-HRP is then introduced, forming a complex with the biotin-labeled antibodies. Then, TMB substrate is added, producing a blue color when it interacts with HRP. After the blue color fully develops, the reaction is terminated with an acidic Stop solution, which changes the blue color to yellow. The intensity of the yellow color is measured at an absorbance of 450 nm and is used to determine the hormone level in the sample.

Materials Required but Not Provided

 Microplate reader capable of measuring absorbance at 450nm.

Materials Provided

- Obesity ELISA Plate (4°C)
- Dilution Buffer (4°C)
- 5x Assay Wash Buffer (4°C)
- Detection Antibody Mix (-20°C)
- Streptavidin-HRP (4°C)
- Substrate (4°C)
- Stop Solution (4°C)

^{**}Spin down small tubes before starting experiment. **

Plasma Sample Preparation

- Centrifuge citrated or EDTA-collected blood at 4°C (1,000 x g for 10 minutes) to separate plasma from erythrocytes. Alternatively, blood collected without anticoagulant can be centrifuged to collect
- 2. Transfer the plasma layer to a new tube without disturbing the buffy layer.
- 3. The plasma may be assayed directly or stored away at -80°C.

Cell Culture Medium Preparation

 Cell culture medium can be assayed directly or stored away at -80°C.

Cell Lysate Preparation

- Wash the cells once with PBS before lysing the cells.
- 3. For a 96-well culture plate, add 40 μL of lysis buffer to each well and incubate at room temperature for 10 minutes.
- 4. Pipette the lysis buffer up and down to detach the cells and transfer the cell lysates into a new tube.
- If necessary, homogenize the cell lysates with a sonicator.
- The cell lysates may be assayed directly or stored at -80°C.
- Use PBS to dilute the cell sample to the appropriate concentration for each assay, if necessary.

Tissue Sample Preparation

- 1. Weigh tissue sample and add 1 mL of lysis buffer per 100mg of tissue.
- 2. Homogenize the tissue samples with a tissue grinder.
- 3. If necessary, further homogenize the tissue samples with a sonicator.
- 4. Centrifuge the sample at 10,000 RPM for 5 minutes to pellet the tissue debris.
- Collect the supernatant and measure the protein concentration of the supernatant. The tissue sample can be assayed directly or stored at -80°C.
- Use PBS to dilute the tissue sample to the appropriate concentration for each assay, if necessary.

Assay Procedure

Reagent preparation before starting experiment

- Dilute the 5x Assay wash buffer to 1x buffer
 - 40 ml 5x Assay wash buffer
 - 160 ml ddH2O.
- Dilute Detection antibody mix 50 times with 1x Diluent buffer.
- Dilute streptavidin-HRP 200 times with 1x Diluent buffer.
- Add 100 μL of sample to each well and incubate for 1 hour at room temperature with gentle shaking.
- Aspirate each well and wash by adding 200 µL of 1x Assay wash buffer. Repeat the process three times for a total of three washes. Make sure all liquid is removed after each wash.
- Add 100 μL of diluted Detection antibody mix to each well and incubate for 1 hour at room temperature with gentle shaking.
- 4. Repeat the aspiration/wash step in step 2.
- Add 100 μL of diluted streptavidin-HRP to each well and incubate for 45 minutes at room temperature with gentle shaking.
- 6. Repeat the aspiration/wash step in step 2.
- 7. Add 100 μL of Substrate to each well and incubate for 10-30 minutes. Substrate incubation time may vary due to differences in antibody reactivity. Stronger blue signals can be stopped earlier. Weaker signals can be incubated for a longer time up to 1 hour. Always stop the reactions of all samples from the same row at the same time.
- 8. Add 50 μ L of Stop solution to each well. The color in the wells should change from blue to yellow.
- 9. Immediately measure the absorbance of the plate at 450 nm using a plate reader.

Obesity Hormone/Adipokine ELISA

	1	2	3	4	5	6	7	8	9	10	11	12
A	Insulin											
В	Amylin											
С	Glucagon											
D	GLP-1											
Е	Leptin											
F	Adiponectin											
G	IL-6											
Н	TNFα											